




Mapping for Moxel Chronic 32 Channel Probe

Moxel connector ASSY-196 – E-1

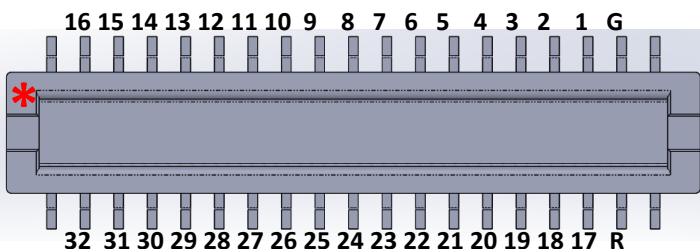


Electrodes are located on the side of the probe as indicated by the arrow below - use Shank A orientation as a reference to construct the channel-map in to your headstage pre-amplifier.

Red Wire and White Wire correspond to indicated REF and GND pins on the probe connector - verify how these connect to your headstage pre-amplifier (often 1 wire is sufficient).

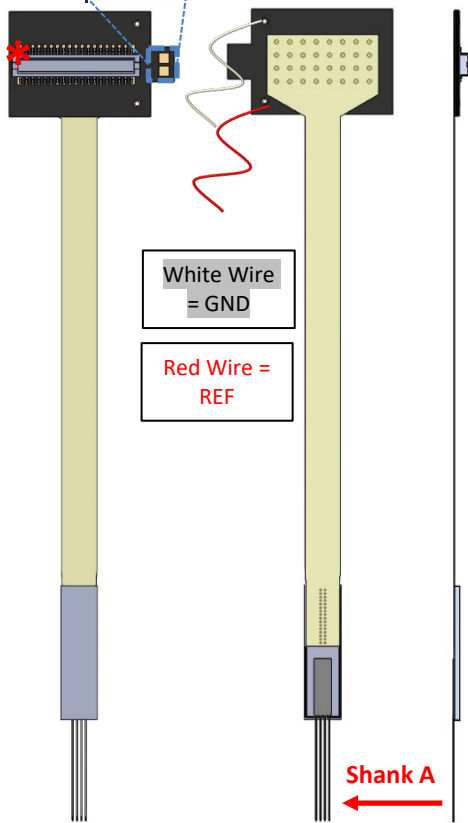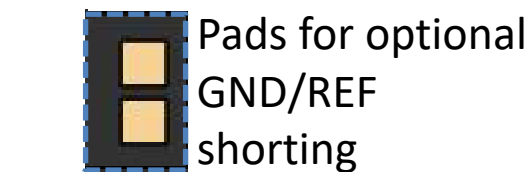

Back view

Front view

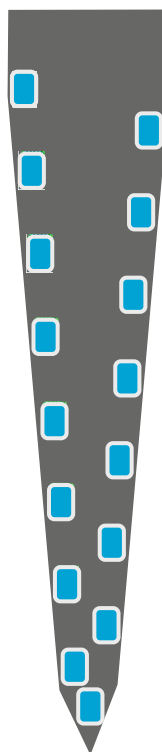

Shank A

31		17	
	32		18
16		2	
	15		1
29		19	
	30		20
14		4	
	13		3
27		21	
	28		22
12		6	
	11		5
9		7	
	25		23
16		24	
	10		8