




Mapping for Molex Chronic 32 Channel Probe

Molex connector ASSY-196 – P-1

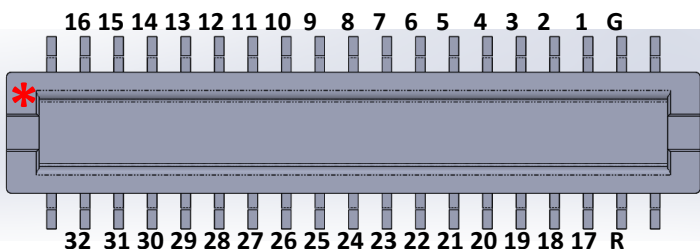

Electrodes are located on the side of the probe as indicated by the arrow below - use Shank A orientation as a reference to construct the channel-map in to your headstage pre-amplifier.


Red Wire and White Wire correspond to indicated REF and GND pins on the probe connector - verify how these connect to your headstage pre-amplifier (often 1 wire is sufficient).

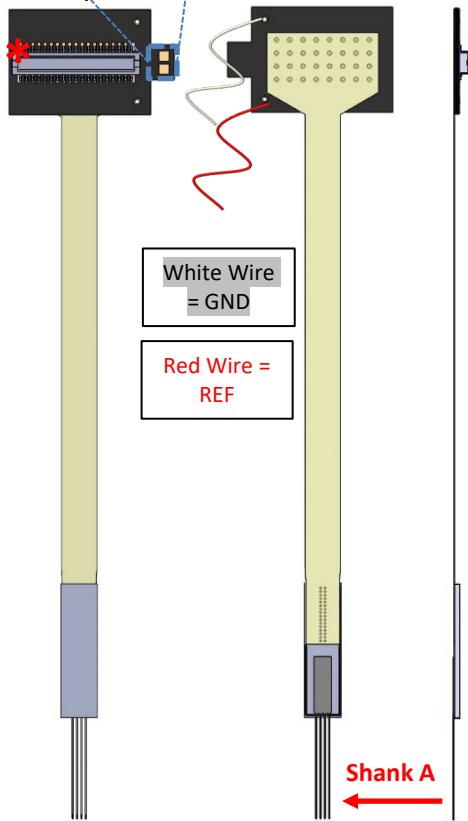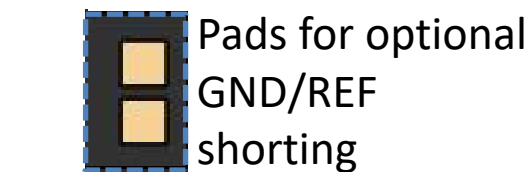

Back view

Front view

Shank A

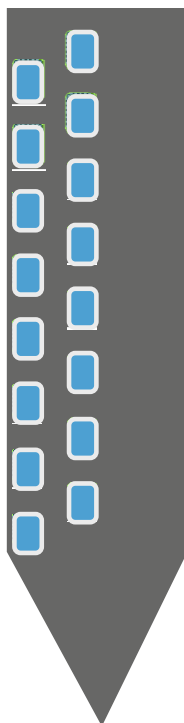

	12		6
31	9	17	7
16	26	2	24
29	10	19	8
14	25	4	23
27	11	21	5
28	13	22	3
30	15	20	1
32		18	