


Mapping for Chronic 64-Channel Probe



ZIF Connector ASSY-236 M2 (June 2022 onwards)

Red Wire and White Wire correspond to indicated REF and GND pins on the probe connector - verify how these connect to your headstage pre-amplifier (1 wire may be sufficient if you prefer GND / REF shorted-together in your headstage).

TOP

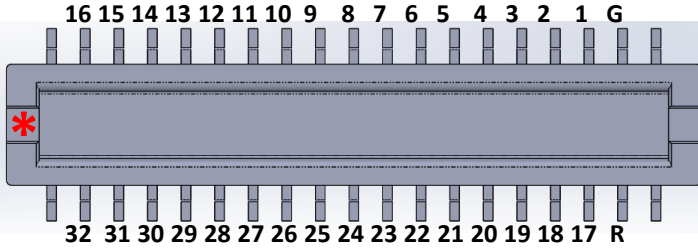

BOTTOM

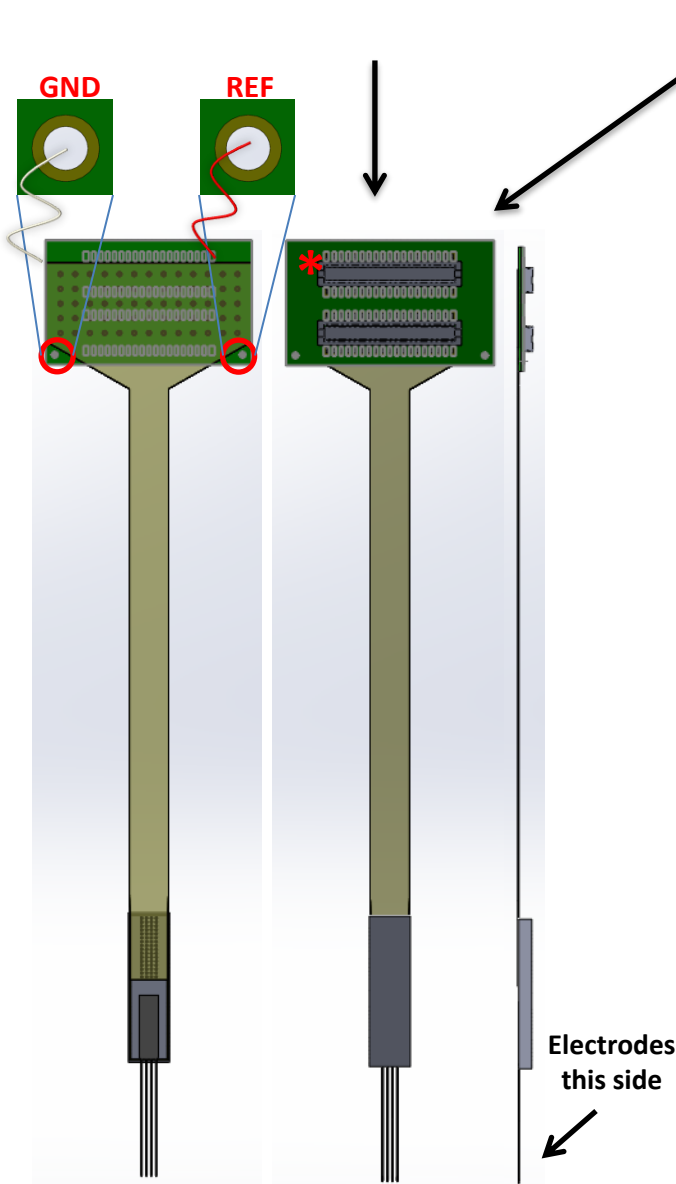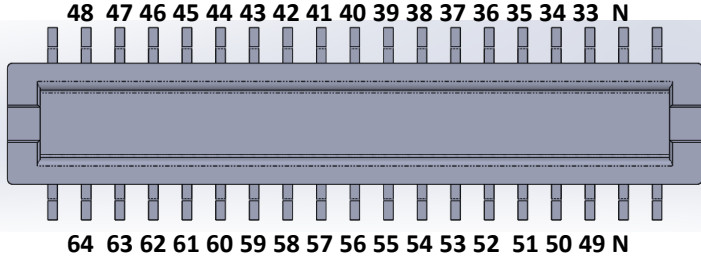

64	34
48	18
16	54
61	23
13	7
60	38
29	22
12	40
59	56
32	8
47	41
14	55
31	20
62	36
15	6
63	39
46	49
30	19
58	1
27	35
10	4
42	52
26	21
24	5
25	53
9	17
43	33
57	3
28	37
44	51
11	2
45	50

Upper-half

Lower-half

